



verity[®] 4060 and 4260

Single and Dual Syringe Pumps

User's Guide



Gilson, Inc. | PO Box 620027 | Middleton, WI 53562-0027 | Tel: 608-836-1551 OR 800-445-7661 | Fax: 608-831-4451

Gilson S.A.S. | 19, avenue des Entrepreneurs | BP 145, F-95400 Villiers-le-bel, FRANCE

www.gilson.com | sales@gilson.com | service@gilson.com | training@gilson.com

LT801540-01

Table of Contents

Safety

Electronic and Hazard Symbols.....	2
Safety Notices	2
Voltage.....	3
Stacking Bracket.....	3
Solvents.....	3
Replacement Parts.....	3

Sécurité

Symboles Électroniques et de Dangers	6
Notes de Sécurité.....	6
Tension.....	7
Kit de superposition	7
Solvants.....	7
Pièces détachées	7

1 Introduction

Description.....	10
Unpacking	11
Standard Equipment.....	11
Accessories.....	12
Customer Service.....	12
Trademarks.....	12
Technical Specifications.....	13

2 Installation

Syringe Installation	16
250 μ L and 500 μ L Syringes	16
All Syringes.....	16
Plumbing Connections	17
GX-241 ASPEC™ or GX-271 ASPEC™ with VERITY® 4060 Single Syringe Pump	17
GX-271 ASPEC™ with VERITY® 4260 Dual Syringe Pump	25
GX-274 ASPEC™ with two VERITY® 4260 Dual Syringe Pumps	32
Rear Panel Connections.....	40
Rear Panel Diagram.....	40
USB Device Port	40
Power Connection	40
Stacking Bracket Installation	41

3 Operation

Front Panel	44
Power Indicator Light	44
Error Indicator Light	44
Side Panel	44
Power Switch	44
Start Up	45
Prime the Syringe Pump	45

4 Maintenance

Helpful Hints	48
Cleaning	49
Exterior.....	49
Syringe.....	49
Valve	51
Part Replacement.....	52
Tubing	52
Syringe.....	52
Valve	53

5 Troubleshooting

Mechanical	56
Electrical	56
Hydraulic.....	56
Error Codes and Messages	58
Repair and Return Policies	59
Before Calling Us.....	59
Warranty Repair	59
Non-Warranty Repair.....	59
Return Procedure	59
Unit End-of-Life.....	59

A Parts and Accessories

GX-Series ASPEC Systems	61
GX-241 ASPEC™ with VERITY® 4060 Single Syringe Pump	61
GX-271 ASPEC™ with VERITY® 4060 Single Syringe Pump	61
GX-271 ASPEC™ with VERITY® 4260 Dual Syringe Pump	61
GX-274 ASPEC™ with Two VERITY® 4260 Dual Syringe Pumps	61
Syringes	62
Piston Seals	62
Valve and Valve Parts	62
Tubing	63
Inlet	63
Vent	63
Plumbing Packages	63
SPE Gas Pressure Regulator	63
Cables and Power Cords	64

B Materials

Liquid Contact Materials	65
--------------------------------	----

Safety

Read this chapter before installing and operating the instrument.

Only trained technical personnel in a laboratory environment may use the instrument for non-medical, liquid handling purposes. For safe and correct use of the instrument, operating and service personnel must follow all instructions contained in this guide when installing, cleaning, and maintaining the instrument. All safety precautions must be observed during all phases of operation, service, and repair of the instrument.

Failure to comply with these precautions or with warnings described in the user's guide violates safety standards of design, manufacture, and intended use of the instrument. Gilson assumes no liability for customers failing to comply with these requirements.

The instrument has been certified to safety standards required in Canada, Europe, and the United States. Refer to the rear panel label on the instrument and the Declaration of Conformity document for the current standards to which the instrument has been found compliant.

Electronic and Hazard Symbols

The following electronic and hazard symbols may appear on the instrument:

Symbol	Explanation
	Direct current Courant continu Gleichstrom
	Protective conductor terminal Borne de terre de protection Schutzleiteranschluss
	Electrical power ON Sous tension Netzschalter ein
	Electrical power OFF Hors tension Netzschalter aus
	Caution Attention Vorsicht
	Caution, risk of electric shock Attention, risque de choc électrique Vorsicht, Elektroschockgefahr
	Caution, hot surface Attention, surface chaude Vorsicht, heiße Oberfläche

Safety Notices

The following safety notices may appear in this document:

 WARNING	WARNING indicates a potentially hazardous situation which, if not avoided, may result in serious injury
 CAUTION	CAUTION indicates a potentially hazardous situation which, if not avoided, may result in minor or moderate injury
 NOTICE	NOTICE indicates a potentially hazardous situation which, if not avoided, may result in equipment damage

Voltage

Ensure that the rear panel is easily accessible. Detach all sources of voltage from the instrument before the service, repair, or exchange of parts. Use only the grounded AC cord provided. Ungrounded power cords can result in electrical shock and serious personal injury. Use only approved fuses with the specified current rating. The instrument must be operated within the voltage specified on the rear panel of the instrument.

Stacking Bracket

The supplied stacking bracket must be used to prevent tipping when stacking two VERITY® 4260 Dual Syringe Pumps. Do not stack more than one pump on top of another pump. The stacking bracket stabilizes the pump stack by anchoring it to the GX-274 ASPEC™.

Solvents

Observe safe laboratory practices when handling solvents. Adequate safety precautions, such as proper ventilation, safety glasses, etc., must be used when handling dangerous liquids. Refer to the Material Safety Data Sheet (MSDS) for each solvent before use.

Replacement Parts

Only use the replacement parts mentioned in this user's guide.

Sécurité

Merci de lire attentivement cette section avant toute installation ou utilisation de l'instrument.

Cet instrument est exclusivement destiné à être utilisé dans un environnement de laboratoire, par un personnel qualifié, à des fins de manipulations de liquides non-médicales. Pour une utilisation correcte et en toute sécurité de l'instrument, il est nécessaire que le personnel qui utilise et réalise la maintenance de l'instrument, suive les instructions contenues dans ce guide lors de l'installation, du nettoyage et de la maintenance de l'instrument. Toutes les consignes de sécurité doivent être respectées durant toutes les phases de fonctionnement, d'entretien ou de réparation de l'instrument.

Le non-respect de ces précautions ou des avertissements spécifiques mentionnés dans ce guide compromet les normes de sécurité de conception, de fabrication et d'utilisation prévue de l'instrument. Gilson décline toute responsabilité en cas d'incapacité du client à se conformer à ces exigences.

L'instrument a été certifié conformément aux normes de sécurité en vigueur au Canada, en Europe et aux Etats-Unis. Merci de vous reporter aux indications mentionnées sur le panneau arrière de l'instrument ainsi qu'au document de Déclaration de Conformité aux normes pour lesquelles l'instrument a été déclaré conforme.

Symboles Électroniques et de Dangers

Les symboles électroniques et de dangers suivants peuvent apparaître sur l'instrument:

Symbol	Explanation
	Direct current Courant continu Gleichstrom
	Protective conductor terminal Borne de terre de protection Schutzleiteranschluss
	Electrical power ON Sous tension Netzschalter ein
	Electrical power OFF Hors tension Netzschalter aus
	Caution Attention Vorsicht
	Caution, risk of electric shock Attention, risque de choc électrique Vorsicht, Elektroschockgefahr
	Caution, hot surface Attention, surface chaude Vorsicht, heiße Oberfläche

Notes de Sécurité

Les notes de sécurité suivantes peuvent apparaître dans ce document:

 WARNING	WARNING indique une situation potentiellement dangereuse qui, si elle n'est pas évitée, peut entraîner des blessures graves
 CAUTION	CAUTION indique une situation potentiellement dangereuse qui, si elle n'est pas évitée, peut entraîner des blessures mineures à modérées
 NOTICE	NOTICE indique une situation potentiellement dangereuse qui, si elle n'est pas évitée, peut entraîner des dégâts matériels

Tension

S'assurer que l'accès au panneau arrière est libre. Déconnecter la source d'alimentation avant toute opération d'entretien, de réparation ou de remplacement de pièces. Utiliser exclusivement le bloc et le cordon d'alimentation avec raccordement à la terre fournis. Un cordon d'alimentation sans terre peut provoquer choc électrique et graves blessures. Utiliser exclusivement des fusibles de l'intensité et du type spécifié. Pour le fonctionnement, respecter la tension indiquée sur le panneau arrière de l'instrument.

Kit de superposition

Le kit de superposition fourni doit être utilisé afin d'éviter tout risque de basculement lors de l'empilage de deux pompes bi-seringues VERITY® 4260. Ne pas empiler plus de 2 pompes. Le kit de superposition doit être fixé au GX-274 ASPEC™.

Solvants

Respecter les Bonnes Pratiques de Laboratoire lors de la manipulation de solvants. Si des liquides dangereux sont utilisés, s'assurer que la ventilation est adéquate et porter en permanence un équipement de protection individuelle (EPI), tel que : lunettes, gants et vêtements de protection. Se reporter aux Fiches de Données de Sécurité relatives aux solvants avant toute utilisation.

Pièces détachées

Utiliser exclusivement les pièces détachées mentionnées dans le guide de l'utilisateur.

Introduction

Chapter One

This chapter provides information on the following topics:

- **Description** on page 10
- **Unpacking** on page 11
- **Customer Service** on page 12
- **Trademarks** on page 12
- **Technical Specifications** on page 13

Description

The VERITY® 4060 Single Syringe Pump and VERITY® 4260 Dual Syringe Pump when paired with a GX-Series ASPEC™, can automate liquid handling procedures and solid phase extraction (SPE) using positive pressure elution in cartridge-based applications.

The VERITY® 4060 Single Syringe Pump is equipped with a user-selectable, small- or large-capacity syringe, a valve for directing liquid from reservoir or probe or air from an external gas source, and an independent pressure-sensing module that monitors pressure on the fluid path that can be used to handle error conditions, such as high cartridge pressure or clogged liquid lines. An SPE air push can be performed on SPE cartridges using the syringe or an external gas source.

The VERITY® 4260 Dual Syringe Pump is equipped with up to two syringes that are user-selectable, can be the same or different, and can be a small- or large-capacity. Two valves direct liquid from reservoirs or up to two probes, or direct air from an external gas source. Two independent pressure-sensing modules monitor pressure on the fluid paths and can be used to handle error conditions, such as high cartridge pressure or clogged liquid lines. An SPE air push can be performed on SPE cartridges using the syringes or an external gas source. When two VERITY® 4260 Dual Syringe Pumps are used with a GX-274 ASPEC™, up to four fluid paths can be controlled and monitored independently, which allows for a high throughput configuration capable of processing up to four samples in parallel.

The following configurations are supported:

- GX-241 ASPEC™ System with VERITY® 4060 Single Syringe Pump
- GX-271 ASPEC™ System with VERITY® 4060 Single Syringe Pump
- GX-271 ASPEC™ System with VERITY® 4260 Dual Syringe Pump
- GX-274 ASPEC™ System with two VERITY® 4260 Dual Syringe Pumps



Figure 1: GX-274 ASPEC™ with Two VERITY® 4260 Dual Syringe Pumps

Unpacking

The syringe pump(s) are delivered with most major components already assembled. Keep the original container and packing assembly in case the syringe pump(s) must be returned to the factory.

To unpack each syringe pump:

1. Open the box and then remove the envelope inside.
2. Lift the syringe pump and foam inserts out of the box and then remove the foam inserts.
3. Remove the cardboard insert containing the accessories and the power supply.

Standard Equipment

Once the syringe pump and the accessories have been unpacked, you should have the following:

- VERITY® 4060 Single Syringe Pump or one or two VERITY® 4260 Dual Syringe Pumps
- Accessory kit that includes:
 - Power cords
 - USB cable
 - Valve key
 - Vent tubing
- Power supply

Documentation

The following documents are included with the syringe pump:

- Declaration of Conformity
- Hazardous Materials Declaration (China RoHS)
- Installation Qualification (IQ) procedure
- Items Included Checklist
- Quality Control (QC) checklist
- Setup Overview

This user's guide (and other ASPEC™ System user's guides), plumbing diagrams, and IQ procedure documents are provided on the ASPEC™ Systems Documentation CD provided with the GX-series ASPEC.

Accessories

Required

Some accessories are required, but are ordered separately:

- Plumbing package
- Syringes

Optional

The following optional accessories are available:

- Air/gas plumbing package
- SPE gas pressure regulator

Customer Service

Gilson, Inc. and its worldwide network of representatives provide customers with the following types of assistance: sales, technical support, applications, and instrument repair.

If you need assistance, please contact your local Gilson representative. Specific contact information can be found at www.gilson.com. To help us serve you quickly and efficiently, please refer to **Before Calling Us** on page 59.

Trademarks

The following trademark may appear in this document:

- Ekonol® is a registered trademark of Saint Gobain Advanced Ceramics LLC.
- Teflon® is a registered trademark of E.I. du Pont de Nemours & Co, Inc.

Technical Specifications

Please be aware of the following before operating the instrument.

NOTICE

Changes or modifications to the instrument not expressly approved by Gilson could void the factory-authorized warranty.

This instrument complies with part 15 of the FCC Rules. Operation is subject to the following two conditions: (1) this instrument may not cause harmful interference, and (2) this instrument must accept any interference received, including interference that may cause undesired operation.

Shielded cables must be used with the instrument to ensure compliance with the FCC Class A limits.

VERITY® 4060 Single and VERITY® 4260 Dual Syringe Pumps

Technical Specification	Definition	
Communication	USB	
Dimensions (W x D x H)	VERITY® 4060 Single Syringe Pump 14.6 x 17.1 x 29.5 cm (5.8 x 6.7 x 11.6 in.) VERITY® 4260 Dual Syringe Pump 22.6 x 17.1 x 29.5 cm (8.9 x 6.7 x 11.6 in.)	
Environmental Conditions	Indoor use Altitude: up to 2000 m Temperature range: 5°C–40°C Humidity: Maximum relative humidity 80% for temperatures up to 31° C, decreasing linearly to 50% relative humidity at 40°C	
Front Panel	Indicator lights for power and error	
Liquid Contact Materials* <small>*For more information, refer to the Materials appendix.</small>	<i>Description</i>	<i>Material</i>
	Syringe	Glass PTFE Ekonol
	Syringe Valve	PEEK Ceramic PTFE
<i>VERITY® 4060 Single and VERITY® 4260 Dual Syringe Pumps Technical Specifications (continued on page 14)</i>		

VERITY® 4060 Single and VERITY® 4260 Dual Syringe Pumps

Maximum Syringe Flow Rate for Water	Syringe Size	Recommended Maximum Flow Rate
	250 µL	10 mL/min
	500 µL	20 mL/min
	1 mL	40 mL/min
	5 mL	100 mL/min
	10 mL	100 mL/min
	25 mL	100 mL/min
Power Requirements	<p>VERITY® 4060 Single Syringe Pump Voltage: 24V DC Current Rating: 1A, 24W</p> <p>VERITY® 4260 Dual Syringe Pump Voltage: 24V DC Current Rating: 2A, 48W</p> <p>External Power Supply Voltage Input Frequency: 50 to 60 Hz Voltage: 100–240V AC Voltage Output Voltage: 24V DC Current Rating: 2.5A, 60W</p>	
Safety and Compliance	<p>The instrument has been certified to safety standards specified for Canada, Europe, and the United States. Refer to the instrument rear panel label and the Declaration of Conformity document for the current standards to which the instrument has been found compliant.</p>	
Syringe Capacity	250 µL, 500 µL, 1 mL, 5 mL, 10 mL, or 25 mL	
Software Control	PC control via USB and TRILUTION® LH Software	
Volumetric Accuracy*	±2% (10%–90% syringe capacity, water)	
Weight	<p>VERITY® 4060 Single Syringe Pump 4.7 kg (10.4 lbs.)</p> <p>VERITY® 4260 Dual Syringe Pump 7.6 kg (16.7 lbs.)</p>	
<p><i>VERITY® 4060 Single and VERITY® 4260 Dual Syringe Pumps Technical Specifications (continued on page 14)</i></p>		

Installation

Chapter Two

This chapter provides information on the following topics:

- [Syringe Installation](#) on page 16
- [Plumbing Connections](#) on page 17
- [Rear Panel Connections](#) on page 40
- [Stacking Bracket Installation](#) on page 41

Syringe Installation

CAUTION To prevent injury, when operating the syringe pump, keep hands clear of syringe to avoid risk of personal injury by pinching.

250 μ L and 500 μ L Syringes

The 250 μ L and 500 μ L syringes are supplied with a cover seal to ensure an airtight fit between the syringe and the valve.

Before mounting one of the smaller syringes (250 μ L, 500 μ L and 1 mL), manually prime it. Place its open end in the pump reservoir and use the piston to aspirate the liquid. This manual prime is not necessary for the 5 mL, 10 mL, and 25 mL syringes.

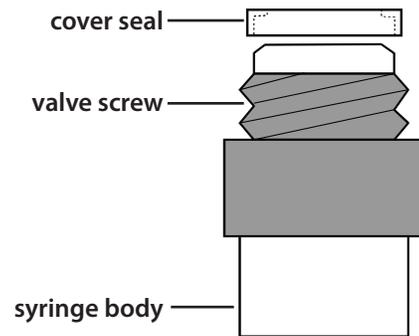


Figure 2: Cover Seal on Syringe

All Syringes

The piston operating rod will be shipped in the down position. If the rod is not in the down position, refer to the instructions for replacing a syringe on page 21. Those instructions include how to lower the rod.

The following procedure is important for correct syringe piston alignment. Improper alignment may cause premature piston failure.

1. Remove the valve from the front panel by removing the two securing valve screws.
2. Lubricate the piston with an alcohol-based solvent (methanol, for example) to reduce piston seal friction during syringe installation.
3. Remove the cap covering port 0 and then loosely screw the syringe into the valve. Do not fully tighten.
4. Loosely attach the valve to the syringe pump with the supplied screws.
5. Pull down the piston so it comes into contact with the piston operating rod, and firmly tighten the piston holding screw.
6. Fully tighten the valve screws to secure the valve.
7. Fully tighten the syringe to the valve.

Remember or note the size of the syringe installed for later software configuration.

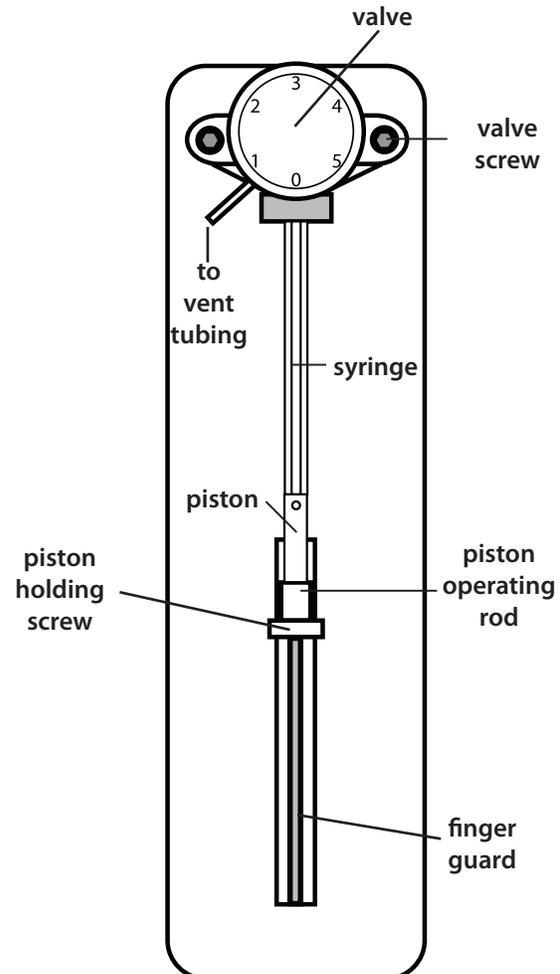


Figure 3: Syringe Installation

Plumbing Connections

GX-241 ASPEC™ or GX-271 ASPEC™ with VERITY® 4060 Single Syringe Pump

Before making the plumbing connections, locate the plumbing package (ordered separately). The plumbing package contains:

- Solvent inlet tubing
- Transfer tubing
- Pressure sensor to valve tubing assembly
- Pump to probe tubing assembly

If the air push will be performed with air/gas (not syringe), locate the air/gas plumbing package (ordered separately, part number 2644703) that contains tubing and fittings. Instructions for [Air/Gas Configuration](#) begin on page 18.

Instructions for [No Air/Gas Configuration](#) begin on page 22.

For instructions to make plumbing connections for [GX-271 ASPEC™ with VERITY® 4260 Dual Syringe Pump](#), refer to the instructions beginning on page 25.

For instructions to make plumbing connections for [GX-274 ASPEC™ with two VERITY® 4260 Dual Syringe Pumps](#), refer to the instructions beginning on page 32.

Air/Gas Configuration

Refer to the diagram and table for instructions to make the plumbing connections for an air/gas configuration.

Plumbing Connections for GX-241 ASPEC™ or GX-271 ASPEC™ with VERITY® 4060 Single Syringe Pump Air/Gas Configuration (with SPE Gas Pressure Regulator)			
Valve	Tubing	Connection	Diagram
Port 1 to pressure sensor OUT	Pressure sensor to valve tubing (part number 4996752)		
	4.5 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	Connect one end to port 1 on the syringe pump and the other end to pressure sensor OUT.	
Port 2 from reservoir	Solvent inlet tubing (part number 499484021)		
	40 inches of Teflon® tubing (0.085" ID x 1/8" OD)	Connect the end of the tubing with the fitting attached to port 2 on the syringe pump. Place the other end in a reservoir.	
Port 3 to gas	GX-241/GX-271 ASPEC Air/Gas Plumbing Package (part number 2644703)		
	Polyurethane tubing (5/64" ID x 1/8" OD)	<p>Install the nut and ferrule to one end of the tubing and connect it to port 3 on the syringe pump. Refer to the diagram on page 21.</p> <p>Connect the other end of the tubing to the SPE gas pressure regulator (part number 25051376). Refer to Air/Gas Supply Connections on page 21.</p>	

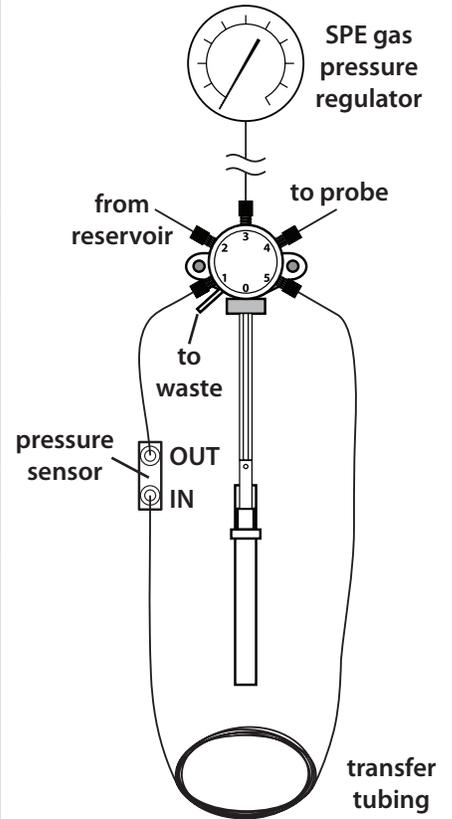


Figure 4: Plumbing Connections for Air/Gas Configuration (with SPE Gas Pressure Regulator)

Plumbing Connections for GX-241 ASPEC™ or GX-271 ASPEC™ with VERITY® 4060 Single Syringe Pump Air/Gas Configuration (with SPE Gas Pressure Regulator)

Valve	Tubing	Connection	Diagram
Vent to waste	Vent tubing (part number F4420577)		
	FEP tubing	Connect one end directly to the vent and place the other end in a suitable receptacle. The valve releases liquid through the waste outlet if the pressure inside the valve is too high.	
Port 4 to probe	Pump to probe tubing (part number 49967722)		
	72 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	Connect the end with the nut and ferrule to port 4 and connect the end with the headless nut to the probe.	

Plumbing Connections for GX-241 ASPEC™ or GX-271 ASPEC™ with VERITY® 4060 Single Syringe Pump Air/Gas Configuration (with SPE Gas Pressure Regulator)

Valve	Tubing	Connection	Diagram
Port 5 to pressure sensor IN	<p>Note: It is recommended to store the excess transfer tubing in the compartment under the syringe pump.</p>		
	<p>5 mL transfer tubing (part number 499671142)</p>		
	101 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	Connect one end to port 5 on the syringe pump and the other end to pressure sensor IN.	
	<p>10 mL transfer tubing (part number 499672282)</p>		
	202 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	Connect one end to port 5 on the syringe pump and the other end to pressure sensor IN.	
	<p>25 mL transfer tubing (part number 4996757021)</p>		
	505 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	Connect one end to port 5 on the syringe pump and the other end to pressure sensor IN.	
<p>For smaller syringe size, use the following transfer tubing:</p> <p>600 µL transfer tubing (part number 49942607)</p> <p>1.1 mL transfer tubing (part number 499424013)</p>			

Air/Gas Supply Connections

The following steps describe how to connect the air/gas tubing.

The tubing shown is blue polyurethane (5/64" ID x 1/8" OD). Fittings are Upchurch P-331 nut (part number 490410332) and Upchurch P-359 ferrule (part number 49041027).

1. Install the nut and ferrule on one end of the tubing.
2. Connect the other end to the pressure regulator. Refer to [Figure 5: Air/Gas Supply Connections](#) for more information.

INSTALLING THE SPE GAS PRESSURE REGULATOR

The SPE gas pressure regulator (part number 25051376, ordered separately) includes the following:

- Double cable clamp
- Hex nut
- Lock washer
- Pressure regulator with attached bracket
- Truss head screw

To install the pressure regulator:

1. Attach the double cable clamp to the bracket using the screw, hex nut, and washer.
2. Attach the double cable clamp to the cable support rod on the GX-241 ASPEC™ or GX-271 ASPEC™.
3. Use the appropriate tubing and fittings to connect the gas pressure regulator to your gas supply.

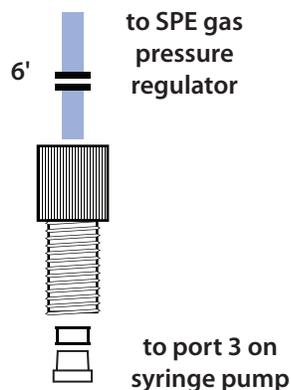


Figure 5: Air/Gas Supply Connections

No Air/Gas Configuration

Refer to the diagram and table for instructions to make the plumbing connections for a no air/gas configuration.

Note: Ports 3, 4, and 5 are not used when there is no air/gas connected. The syringe pump is shipped with a plug in each port. Leave the plug installed in all unused ports.

Plumbing Connections for GX-241 ASPEC™ or GX-271 ASPEC™ with VERITY® 4060 Single Syringe Pump No Air/Gas Configuration			
Valve	Tubing	Connection	Diagram
Port 1 to pressure sensor OUT	<p>Pressure sensor to valve tubing (part number 4996752)</p> <p>4.5 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)</p>	Connect one end to port 1 on the syringe pump and the other end to pressure sensor OUT.	<p>Figure 6: Plumbing Connections for No Air/Gas Configuration</p>

**Plumbing Connections for GX-241 ASPEC™ or GX-271 ASPEC™ with VERITY® 4060 Single Syringe Pump
No Air/Gas Configuration**

Valve	Tubing	Connection	Diagram
Pressure sensor IN to probe	<p>Note: The transfer tubing is coiled. To connect one end of the tubing to the probe, the clips that hold the tubing will need to be cut. After the connection to the probe has been made, it is recommended to re-coil the transfer tubing and place the excess tubing in the compartment under the syringe pump.</p>		
	<p>5 mL transfer tubing (part number 499671142)</p>		
	101 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	Connect one end to pressure sensor IN and the other end to the probe.	
	<p>10 mL transfer tubing (part number 499672282)</p>		
	202 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	Connect one end to pressure sensor IN and the other end to the probe	
	<p>25 mL transfer tubing (part number 4996757021)</p>		
	505 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	Connect one end to pressure sensor IN and the other end to the probe	
<p>For smaller syringe size, use the following transfer tubing:</p> <p>600 µL transfer tubing (part number 49942607)</p> <p>1.1 mL transfer tubing (part number 499424013)</p>			

**Plumbing Connections for GX-241 ASPEC™ or GX-271 ASPEC™ with VERITY® 4060 Single Syringe Pump
No Air/Gas Configuration**

Valve	Tubing	Connection	Diagram
Port 2 from reservoir	<p>Solvent inlet tubing (part number 499484021)</p> <p>40 inches of Teflon® tubing (0.085" ID x 1/8" OD)</p>	<p>Connect the end of the tubing with the fitting attached to port 2 on the syringe pump. Place the other end in a reservoir.</p>	
Vent to waste	<p>Vent tubing (part number F4420577)</p> <p>FEP tubing</p>	<p>Connect one end directly to the vent and place the other end in a suitable receptacle. The valve releases liquid through the waste outlet if the pressure inside the valve is too high.</p>	

GX-271 ASPEC™ with VERITY® 4260 Dual Syringe Pump

Before making the plumbing connections, locate the plumbing package (ordered separately).

The plumbing package contains:

- Solvent inlet tubing
- Transfer tubing
- Pressure sensor to valve tubing assembly
- Pump to probe tubing assembly

If the air push will be performed with air/gas (not syringe), locate the air/gas plumbing package (ordered separately) that contains tubing and fittings. Instructions for [Air/Gas Configuration](#) begin on page 26.

Instructions for [No Air/Gas Configuration](#) begin on page 30.

For instructions to make plumbing connections for a [GX-241 ASPEC™](#) or [GX-271 ASPEC™ with VERITY® 4060 Single Syringe Pump](#), refer to the instructions beginning on page 17.

For instructions to make plumbing connections for a [GX-274 ASPEC™ with two VERITY® 4260 Dual Syringe Pumps](#), refer to the instructions beginning on page 32.

Air/Gas Configuration

Refer to the diagram and table for instructions to make the plumbing connections for an air/gas configuration.

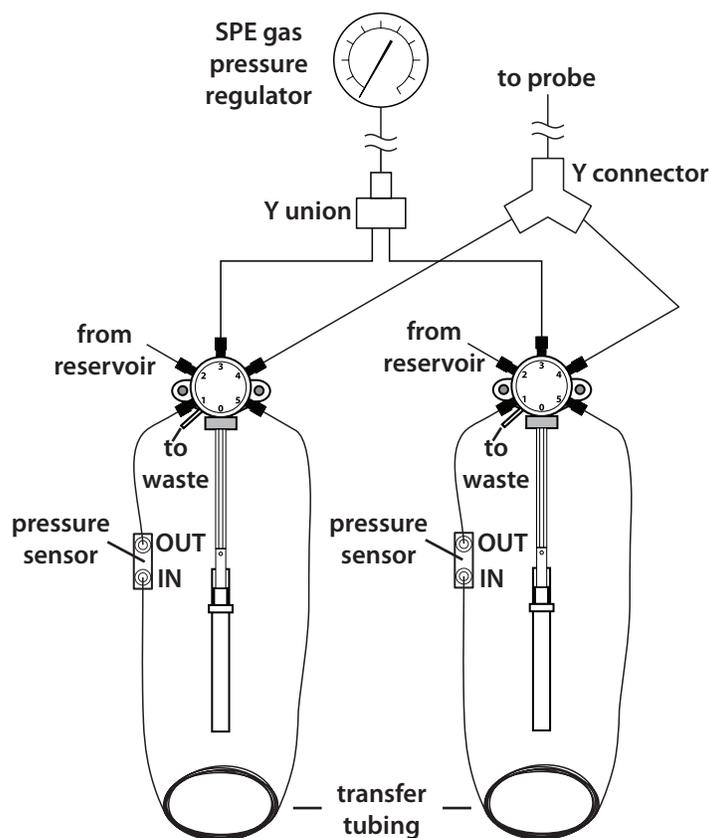


Figure 7: Plumbing Connections for Air/Gas Configuration (with SPE Gas Pressure Regulator)

<i>Plumbing Connections for GX-271 ASPEC™ with VERITY® 4260 Dual Syringe Pump Air/Gas Configuration (with SPE Gas Pressure Regulator)</i>		
<i>Valve</i>	<i>Tubing</i>	<i>Connection</i>
Port 1 to pressure sensor OUT	Pressure sensor to valve tubing (part number 4996752)	
	4.5 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	For each valve, connect one end to port 1 on the syringe pump and the other end to pressure sensor OUT.
Port 2 from reservoir	Solvent inlet tubing (part number 499484021)	
	40 inches of Teflon® tubing (0.085" ID x 1/8" OD)	Connect the end of the tubing with the fitting attached to port 2 on the syringe pump. Place the other end in a reservoir.

Plumbing Connections for GX-271 ASPEC™ with VERITY® 4260 Dual Syringe Pump Air/Gas Configuration (with SPE Gas Pressure Regulator)		
Valve	Tubing	Connection
Port 3 to Y union to gas	GX-271 ASPEC with Dual Syringe Pump Air/Gas Plumbing Package (part number 2644709)	
	Polyurethane tubing (5/64" ID x 1/8" OD)	For each valve, connect the blue polyurethane tubing (with fittings attached) to port 3 on the syringe pump. Connect the remaining end of the blue polyurethane tubing to SPE gas pressure regulator (part number 25051376). Refer to Air/Gas Supply Connections on page 29 for installation instructions.
Port 4 to Y connector	Dual Adapter Kit, GX-271 ASPEC (part number 2644708)	
	24 inches of Teflon® FEP tubing labeled "Dual Adapt" (0.062" ID x 1/8" OD)	There are two lengths of tubing with fittings attached. For each valve, connect one end to port 4 on the syringe pump and connect the other end to the Y connector.
Port 5 to pressure sensor IN	Note: It is recommended to store the excess transfer tubing in the compartments under the syringe pump.	
	5 mL transfer tubing (part number 499671142)	
	101 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	For each valve, connect one end to port 5 on the syringe pump and the other end to pressure sensor IN.
	10 mL transfer tubing (part number 499672282)	
	202 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	For each valve, connect one end to port 5 on the syringe pump and the other end to pressure sensor IN.
	25 mL transfer tubing (part number 4996757021)	
	505 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	For each valve, connect one end to port 5 on the syringe pump and the other end to pressure sensor IN.
For smaller syringe size, use the following transfer tubing: 600 µL transfer tubing (part number 49942607) or 1.1 mL transfer tubing (part number 499424013)		

<i>Plumbing Connections for GX-271 ASPEC™ with VERITY® 4260 Dual Syringe Pump Air/Gas Configuration (with SPE Gas Pressure Regulator)</i>		
<i>Valve</i>	<i>Tubing</i>	<i>Connection</i>
Y connector to probe	Pump to probe tubing (part number 49967722)	
	72 inches of Teflon® FEP tubing (0.062" ID X 1/8" OD)	Connect the end with the nut and ferrule to the Y connector and connect the end with the headless nut to the probe.
Y union to gas	Polyurethane tubing (5/64" ID x 1/8" OD)	Refer to Air/Gas Supply Connections on page 29.
Vent to waste	Vent tubing (part number F4420577)	
	FEP tubing	Connect one end directly to the vent and place the other end in a suitable receptacle. The valve releases liquid through the waste outlet if the pressure inside the valve is too high.

Air/Gas Supply Connections

The following steps describe how to connect the air/gas tubing.

The tubing and Y unions are already assembled.

The tubing shown is blue polyurethane (5/64" ID x 1/8" OD). Fittings are Upchurch P-331 nut (part number 490410332) and Upchurch P-359 ferrule (part number 49041027).

INSTALLING THE SPE GAS PRESSURE REGULATOR

The SPE gas pressure regulator (part number 25051376, ordered separately) includes the following:

- pressure regulator with attached bracket
- truss head screw
- hex nut
- lock washer
- double cable clamp

To install the pressure regulator:

1. Attach the double cable clamp to the bracket using the screw, hex nut, and washer.
2. Attach the double cable clamp to the cable support rod on the GX-271 ASPEC™.
3. Use the appropriate tubing and fittings to connect the gas pressure regulator to your gas supply.

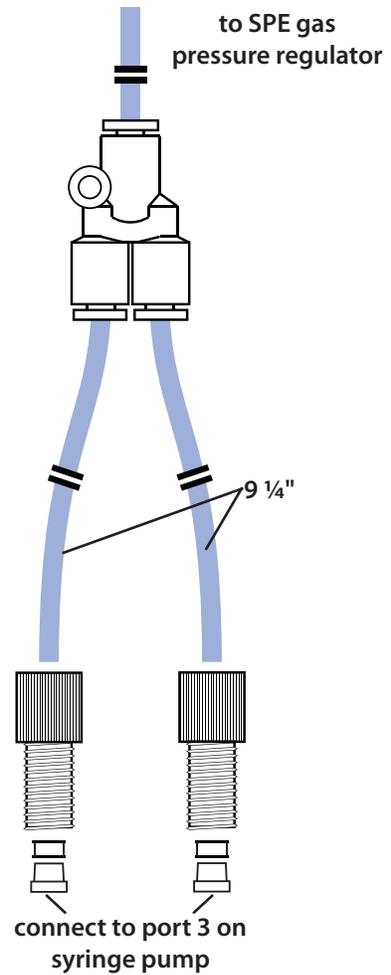


Figure 8: Air/Gas Supply Connections for GX-271 ASPEC™ with VERITY® 4260 Dual Syringe Pump

No Air/Gas Configuration

Refer to the diagram and table for instructions to make the plumbing connections for a no air/gas configuration.

Note: Ports 3, 4, and 5 are not used when there is no air/gas connected. The syringe pump is shipped with a plug in each port. Leave the plug installed in all unused ports.

The Y connector referenced is included in Dual Adapter Kit, GX-271 ASPEC (part number 2644708).

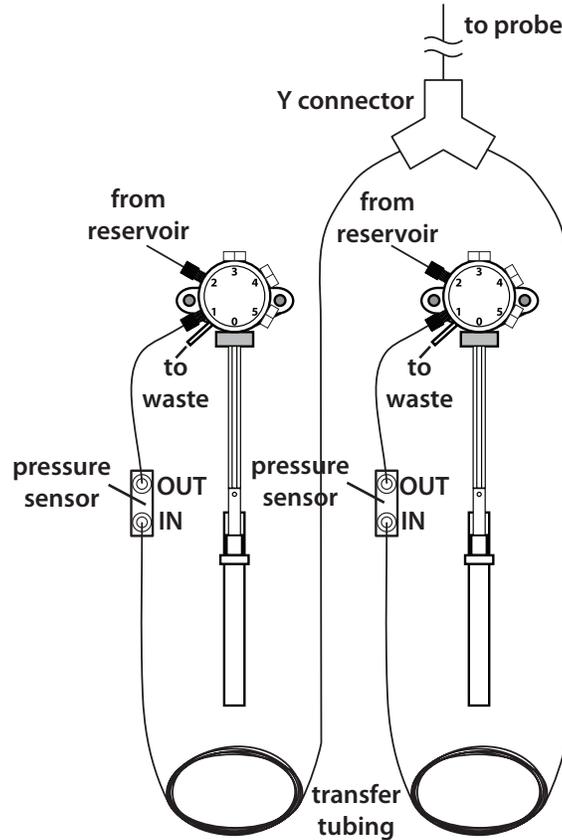


Figure 9: No Air/Gas Configuration

Plumbing Connections for GX-271 ASPEC™ with VERITY® 4260 Dual Syringe Pump No Air/Gas Configuration

Valve	Tubing	Connection
Port 1 to pressure sensor OUT	Pressure sensor to valve tubing (part number 4996752)	
	4.5 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	For each valve, connect one end to port 1 on the syringe pump and the other end to pressure sensor OUT.

Plumbing Connections for GX-271 ASPEC™ with VERITY® 4260 Dual Syringe Pump No Air/Gas Configuration		
Valve	Tubing	Connection
Pressure sensor IN to Y connector	Note: The transfer tubing is coiled. To connect one end of the tubing to the probe, the clips that hold the tubing will need to be cut. After the connection to the probe has been made, it is recommended to re-coil the transfer tubing and place the excess tubing in the compartment under the syringe pump.	
	5 mL transfer tubing (part number 499671142)	
	101 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	For each valve, connect one end to pressure sensor IN and the other end to the Y connector.
	10 mL transfer tubing (part number 499672282)	
	202 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	For each valve, connect one end to pressure sensor IN and the other end to the Y connector.
	25 mL transfer tubing (part number 4996757021)	
	505 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	For each valve, connect one end to pressure sensor IN and the other end to the Y connector.
For smaller syringe size, use the following transfer tubing: 600 µL transfer tubing (part number 49942607) 1.1 mL transfer tubing (part number 499424013)		
Port 2 from reservoir	Solvent inlet tubing (part number 499484021)	
	40 inches of Teflon® tubing (0.085" ID x 1/8" OD)	For each valve, connect the end of the tubing with the fitting attached to port 2 on the syringe pump. Place the other end in a reservoir.
Y connector to probe	Pump to probe tubing (part number 49967722)	
	72 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	Connect the end with the nut and ferrule to the Y connector and connect the end with the headless nut to the probe.
Vent to waste	Vent tubing (part number F4420577)	
	FEP tubing	Connect one end directly to the vent and place the other end in a suitable receptacle. The valve releases liquid through the waste outlet if the pressure inside the valve is too high.

GX-274 ASPEC™ with two VERITY® 4260 Dual Syringe Pumps

Before making the plumbing connections, locate the plumbing package (ordered separately).

The plumbing package contains:

- Solvent inlet tubing
- Transfer tubing
- Pressure sensor to valve tubing assembly
- Pump to probe tubing assembly

If the air push will be performed with air/gas (not syringe), locate the air/gas plumbing package (ordered separately) that contains tubing and fittings. Instructions for [Air/Gas Configuration](#) begin on page 33.

Instructions for [No Air/Gas Configuration](#) begin on page 37.

For instructions to make plumbing connections for [GX-241 ASPEC™](#) or [GX-271 ASPEC™](#) with [VERITY® 4060 Single Syringe Pump](#), refer to the instructions beginning on page 17.

For instructions to make plumbing connections for [GX-271 ASPEC™](#) with [VERITY® 4260 Dual Syringe Pump](#), refer to the instructions beginning on page 25.

Air/Gas Configuration

Refer to the diagram and table for instructions to make the plumbing connections for an air/gas configuration.

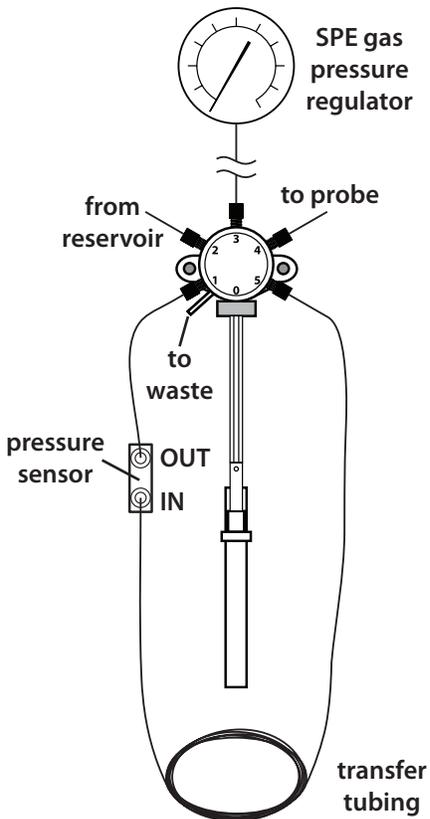
Plumbing Connections for GX-274 ASPEC™ with Two VERITY® 4260 Dual Syringe Pumps Air/Gas Configuration (with SPE Gas Pressure Regulator)			
Valve	Tubing	Connection	Diagram
Port 1 to pressure sensor OUT	Pressure sensor to valve tubing (part number 4996752)		 <p>The diagram shows a syringe pump with a pressure gauge at the top. The gauge is labeled 'SPE gas pressure regulator'. Below the gauge, there are five ports labeled 1 through 5. Port 1 is connected to 'from reservoir'. Port 2 is connected to 'to waste'. Port 3 is connected to 'to probe'. Port 4 is connected to 'pressure sensor OUT'. Port 5 is connected to 'pressure sensor IN'. A 'transfer tubing' is shown at the bottom of the pump assembly.</p>
	4.5 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	For each valve, connect one end to port 1 on the syringe pump and the other end to pressure sensor OUT.	
Port 2 from reservoir	Solvent inlet tubing (part number 499484021)		
	40 inches of Teflon® tubing (0.085" ID x 1/8" OD)	For each valve, connect the end of the tubing with the fitting attached to port 2 on the syringe pump. Place the other end in a reservoir.	

Figure 10: Air/Gas Configuration

**Plumbing Connections for GX-274 ASPEC™ with Two VERITY® 4260 Dual Syringe Pumps
Air/Gas Configuration (with SPE Gas Pressure Regulator)**

Valve	Tubing	Connection	Diagram
Port 3 to gas	GX-274 Air/Gas Plumbing Package (part number 2644707)		
	Polyurethane tubing (5/64" ID x 1/8" OD)	For each valve, connect the blue polyurethane tubing (with fittings attached) to port 3 on the syringe pump. Connect the other end of the tubing to the SPE gas pressure regulator (part number 25051376). Refer to Installing the SPE Gas Pressure Regulator on page 21 for installation instructions.	
Vent to waste	Vent tubing (part number F4420577)		
	FEP tubing	For each valve, connect one end directly to the vent and place the other end in a suitable receptacle. The valve releases liquid through the waste outlet if the pressure inside the valve is too high.	
Port 4 to probe	Pump to probe tubing (part number 49967722)		
	72 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	For each valve, connect the end with the nut and ferrule to port 4 and connect the end with the headless nut to the probe.	

**Plumbing Connections for GX-274 ASPEC™ with Two VERITY® 4260 Dual Syringe Pumps
Air/Gas Configuration (with SPE Gas Pressure Regulator)**

Valve	Tubing	Connection	Diagram
Port 5 to pressure sensor IN	<p>Note: It is recommended to store the excess transfer tubing in the compartment under the syringe pump.</p>		
	<p>5 mL transfer tubing (part number 499671142)</p>		
	101 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	For each valve, connect one end to port 5 on the syringe pump and the other end to pressure sensor IN.	
	<p>10 mL transfer tubing (part number 499672282)</p>		
	202 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	For each valve, connect one end to port 5 on the syringe pump and the other end to pressure sensor IN.	
	<p>25 mL transfer tubing (part number 4996757021)</p>		
	505 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)	For each valve, connect one end to port 5 on the syringe pump and the other end to pressure sensor IN.	
<p>For smaller syringe size, use the following transfer tubing:</p> <p>600 µL transfer tubing (part number 49942607)</p> <p>1.1 mL transfer tubing (part number 499424013)</p>			

Air/Gas Supply Connections

The following steps describe how to connect the air/gas tubing.

The tubing and Y unions are already assembled.

Tubing shown is blue polyurethane (5/64" ID x 1/8" OD). Fittings are Upchurch P-331 nut (part number 490410332) and Upchurch P-359 ferrule (part number 49041027). Tubing lengths indicated assume that the two VERITY® 4260 Dual Syringe Pumps are stacked.

INSTALLING THE SPE GAS PRESSURE REGULATOR

The SPE gas pressure regulator (part number 25051376, ordered separately) includes the following:

- pressure regulator with attached bracket
- truss head screw
- hex nut
- lock washer
- double cable clamp

To install the pressure regulator:

1. Attach the double cable clamp to the bracket using the screw, hex nut, and washer.
2. Attach the double cable clamp to the cable support rod on the GX-274 ASPEC™.
3. Use the appropriate tubing and fittings to connect the gas pressure regulator to your gas supply.

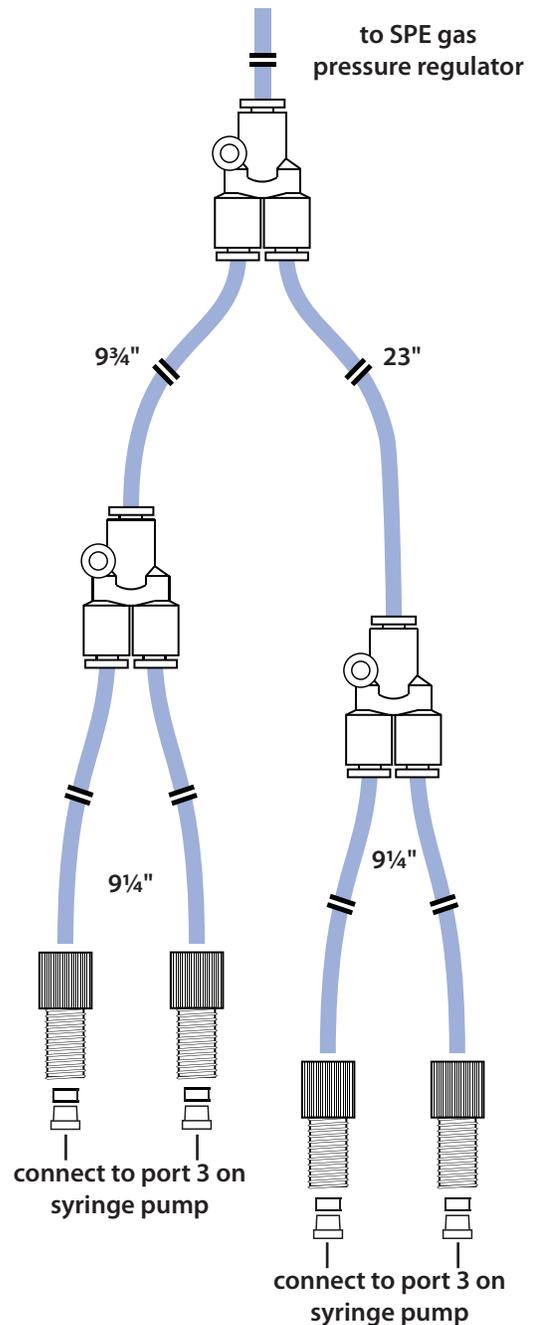


Figure 11: Air/Gas Supply Connections for GX-274 ASPEC™ with Two VERITY® 4260 Dual Syringe Pumps

No Air/Gas Configuration

Refer to the diagram and table for instructions to make the plumbing connections for a no air/gas configuration.

Note: Ports 3, 4, and 5 are not used when there is no air/gas connected. The syringe pump is shipped with a plug in each port. Leave the plug installed in all unused ports.

Plumbing Connections for GX-274 ASPEC™ with Two VERITY® 4260 Dual Syringe Pumps No Air/Gas Configuration			
Valve	Tubing	Connection	Diagram
Port 1 to pressure sensor OUT	<p>Pressure sensor to valve tubing (part number 4996752)</p> <p>4.5 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)</p>	For each valve, connect one end to port 1 on the syringe pump and the other end to pressure sensor OUT.	<p>The diagram shows a syringe pump head with six ports labeled 0 through 5. Port 1 is connected to the 'OUT' port of a pressure sensor. Port 3 is labeled 'from reservoir', port 4 is 'to waste', and port 5 is 'to probe'. Ports 2, 3, 4, and 5 have plugs installed. A pressure sensor is shown with 'IN' and 'OUT' ports. A coil of 'transfer tubing' is connected to the 'to probe' port.</p>

Figure 12: No Air/Gas Configuration

Plumbing Connections for GX-274 ASPEC™ with Two VERITY® 4260 Dual Syringe Pumps
No Air/Gas Configuration

Valve	Tubing	Connection	Diagram
Pressure sensor IN to probe	<p>Note: The transfer tubing is coiled. To connect one end of the tubing to the probe, the clips that hold the tubing will need to be cut. After the connection to the probe has been made, it is recommended to re-coil the transfer tubing and place the excess tubing in the compartment under the syringe pump.</p>		
	<p>5 mL transfer tubing (part number 499671142)</p>		
	<p>101 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)</p>	<p>For each valve, connect one end to pressure sensor IN and the other end to the probe.</p>	
	<p>10 mL transfer tubing (part number 499672282)</p>		
	<p>202 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)</p>	<p>For each valve, connect one end to pressure sensor IN and the other end to the probe</p>	
	<p>25 mL transfer tubing (part number 4996757021)</p>		
	<p>505 inches of Teflon® FEP tubing (0.062" ID x 1/8" OD)</p>	<p>For each valve, connect one end to pressure sensor IN and the other end to the probe</p>	
	<p>For smaller syringe size, use the following transfer tubing:</p> <p>600 µL transfer tubing (part number 49942607)</p> <p>1.1 mL transfer tubing (part number 499424013)</p>		

Plumbing Connections for GX-274 ASPEC™ with Two VERITY® 4260 Dual Syringe Pumps
No Air/Gas Configuration

Valve	Tubing	Connection	Diagram
Port 2 from reservoir	<p>Solvent inlet tubing (part number 499484021)</p> <p>40 inches of Teflon® tubing (0.085" ID x 1/8" OD)</p>	<p>For each valve, connect the end of the tubing with the fitting attached to port 2 on the syringe pump. Place the other end in a reservoir.</p>	<p>The diagram shows a top-down view of a valve assembly. It has six ports labeled 0 through 5. Port 3 is at the top and is labeled 'plug'. Port 4 is to the right and is labeled 'plug'. Port 5 is at the bottom right and is labeled 'plug'. Port 0 is at the bottom and is labeled 'to waste'. Port 1 is at the bottom left. Port 2 is on the left and is labeled 'from reservoir'. A 'pressure sensor' is connected to the valve assembly, with 'OUT' and 'IN' ports. A 'transfer tubing' coil is connected to the 'to waste' port and is labeled 'to probe'. A vertical tube extends downwards from the valve assembly.</p>
Vent to waste	<p>Vent tubing (part number F4420577)</p> <p>FEP tubing</p>	<p>For each valve, connect one end directly to the vent and place the other end in a suitable receptacle. The valve releases liquid through the waste outlet if the pressure inside the valve is too high.</p>	

Rear Panel Connections

Rear Panel Diagram

- 1 USB device port
- 2 Power receptacle

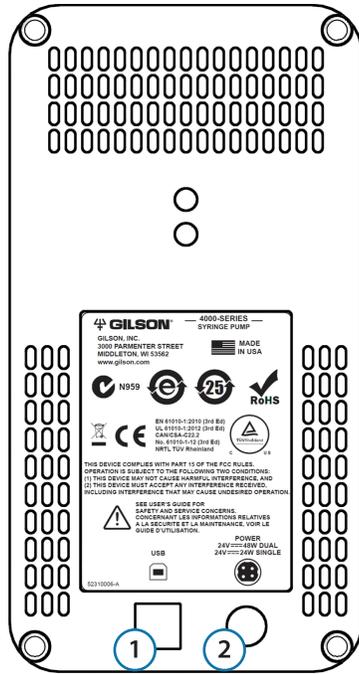


Figure 13: Rear Panel Diagram - VERITY® 4060 Single Syringe Pump

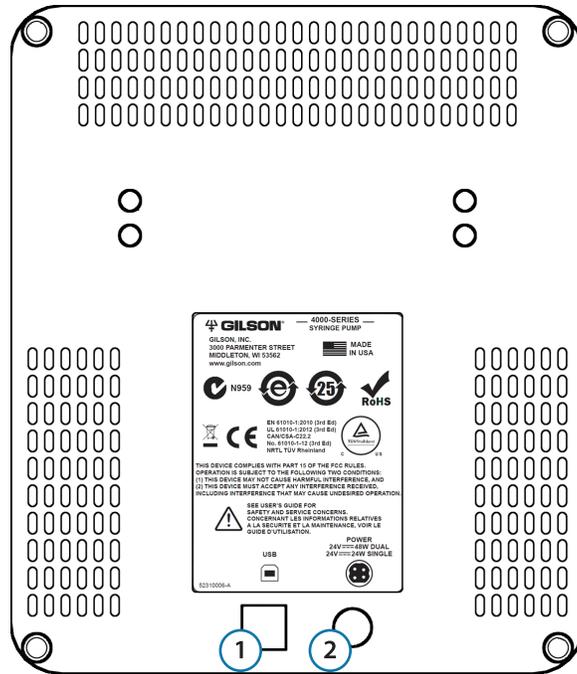


Figure 14: Rear Panel Diagram - VERITY® 4260 Dual Syringe Pump

USB Device Port

The syringe pump communicates with a standard PC via USB.

To make the USB connection between the syringe pump and the controlling device (PC), use the USB cable (part number 38014101) supplied in the accessory kit. Use the end with the “A-type” (flat) connector to connect to the controlling device and use the end with the “B-type” (square) connector to connect to the syringe pump.

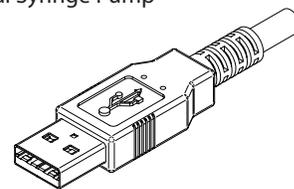


Figure 15: USB Cable with “A-Type” Connector

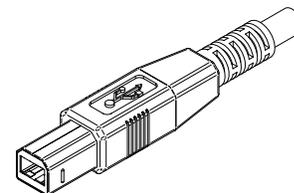


Figure 16: USB Cable with “B-Type” Connector

Power Connection

Use the power cord on the external power supply to make the connection between the power receptacle on the syringe pump and the external power supply.

The connection from the external power supply to the syringe pump uses a connector with a locking collar. Check the alignment of the pins and then push it in until it clicks and locks in place. To disconnect, pull back on the locking collar and then disconnect the cable from the rear panel of the syringe pump.

Locate the appropriate power cord for your line voltage and then make the connection between the external power supply and the AC power source.

Stacking Bracket Installation

The supplied stacking bracket kit (part number 31030017) must be used to prevent tipping when stacking two VERITY® 4260 Dual Syringe Pumps. Do not stack more than one pump on top of another pump. The stacking bracket stabilizes the pump stack by anchoring it to the GX-274 ASPEC™.

The following items will be included with the stacking bracket kit:

- L-bracket
- Two Phillips screws
- Four rubber spacers (pre-installed)
- One metal spacer

To install the stacking bracket:

1. Thread a Phillips screw through the top hole in the L-bracket and through the 0.175" metal spacer.
2. Insert the screw and threaded components into the top hole of the left support on the GX-274 ASPEC. Loosely tighten the top screw.
3. Insert the other screw through the L-bracket into the bottom hole of the left support on the GX-274 ASPEC. Fully tighten the screw.
4. Fully tighten the top screw.
5. Slide one of the VERITY® 4060 Dual Syringe Pump under the installed bracket.
6. Set the second VERITY® 4060 Dual Syringe Pump on top of the first, threading the back support feet through the holes on the L-bracket.



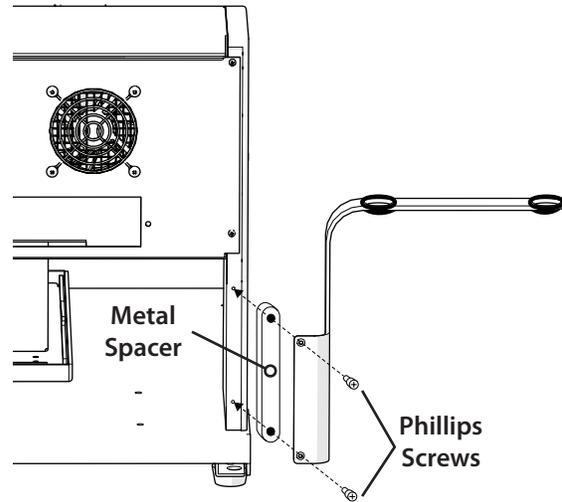



Figure 17: Stacking Bracket and Components (Rear View)

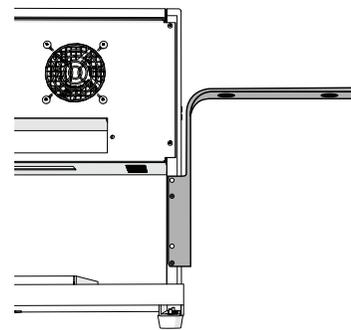


Steps 1 to 4

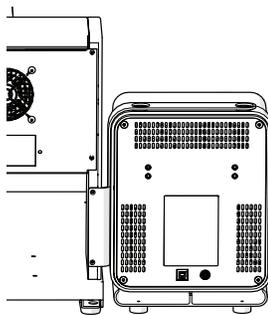

Step 5

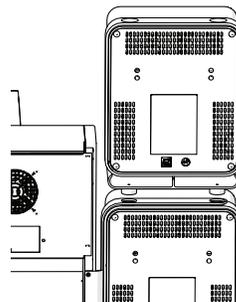

Step 6

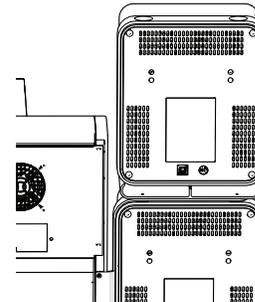

Step 7

Operation

Chapter Three

TRILUTION® LH software provides control of the VERITY® 4060 Single and VERITY® 4260 Dual Syringe Pumps. For more information about TRILUTION LH software, refer to its user's guide and the documentation supplied with the software.

This chapter provides information on the following topics:

- **Front Panel** on page 44
- **Start Up** on page 45
- **Prime the Syringe Pump** on page 45

Front Panel

The front panel includes a power indicator light and an error indicator light.

- ① Power Indicator Light
- ② Error Indicator Light

Power Indicator Light

The green indicator becomes lit when power to the syringe pump is switched on.

Error Indicator Light

The red indicator light flashes when an error has been encountered.

Side Panel

The side panel includes the power switch.

- ③ Power Switch

Power Switch

| indicates that the electrical power is switched on.

O indicates that the electrical power is switched off.



Figure 18: Front/Side Panel Diagram - VERITY® 4060 Single Syringe Pump

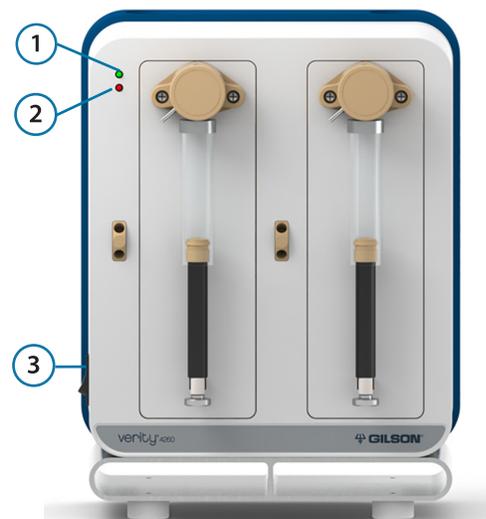


Figure 19: Front/Side Panel Diagram - VERITY® 4260 Dual Syringe Pump

Start Up

To start the syringe pump, turn on power to the syringe pump using the switch on the side panel.

- The power indicator light on the syringe pump front panel illuminates.
- The syringe pump initializes. It stops with the valve set to the outlet (transfer tubing) position.

Prime the Syringe Pump

It is recommended to prime the syringe pump with liquid before using it the first time, or if it has not been used for some time.

The following procedures use TRILUTION® LH Software.

1. Open a method with a VERITY 4060 Syringe Pump or VERITY 4260 Syringe Pump in the configuration.
2. Click **Run**.
3. Add the method to the sample list.
4. Select the Manual Control tab.
5. Select the method from the drop-down list of methods and then click **Go**. The instruments will initialize.
6. If using a VERITY® 4260 Dual Syringe Pump, select the syringe(s) that will be used (Left Syringe and/or Right Syringe).
7. Enter a value for Prime Flow Rate. This is the speed at which reservoir fluid moves into and out of the syringe.

NOTICE

TRILUTION LH does not validate any values or check for safe movement in manual control. Verify entries before clicking any buttons.

8. Click **Prime** to start priming.
9. Click **Stop Prime** to end priming.

Maintenance

Chapter Four

When performing the maintenance described in this chapter, use good laboratory practice, including, but not limited to, wearing protective clothing and preparing the maintenance space for service. After completing the maintenance operation, verify the safe and good working order of the part and instrument.

This chapter contains the following information for maintaining the syringe pump:

- **Helpful Hints** on page 48
- **Cleaning** on page 49
- **Part Replacement** on page 52

Helpful Hints

To keep the syringe pump at peak performance, Gilson recommends the following:

- Change or clean the piston seals and tubing regularly to maintain maximum performance.
- Do not cycle the syringe pump without fluid. Doing this causes excessive piston seal wear.
- Flush the syringe pump daily with distilled or deionized water. On a weekly basis, flush with a 10% solution of bleach or weak detergent.
- If bubbles remain in the syringes after priming, clean the syringes with alcohol.
- Check periodically to ensure that all fittings are tight.
- Check that each syringe is tight in the syringe pump valve.
- Clean the valve if the system has not been used for a while.
- Allow fluids to equilibrate to room temperature before running them through the system; cold fluids may cause leakage.

Cleaning

Exterior

The instrument should be cleaned occasionally using a dry, clean cloth. Or, if necessary, use a cloth dipped in soapy water. If liquid is accidentally spilled on the instrument, wipe it using a dry, clean cloth.

Syringe

Clean a syringe when some or all of the following occurs:

- Corrosive or hazardous liquids have been pumped
- Possible back flow of liquids into the waste tubing
- Leakage
- Aspiration of samples or reagents into the syringe

To clean a syringe, follow the procedures below and use the diagram as a reference.

Remove the Syringe

The following procedures use TRILUTION® LH software.

1. Open a method with a VERITY 4060 Syringe Pump or VERITY 4260 Syringe Pump in the configuration.
2. Click **Run**.
3. Add a method to the sample list.
4. Select the Manual Control tab.
5. Select a method from the drop-down list of methods and then click **Go**. The instruments will initialize.
6. If using a VERITY® 4260 Dual Syringe Pump, select the syringe(s) that will be removed (Left Syringe and/or Right Syringe).
7. Disconnect the syringe piston from the piston operating rod by unscrewing the piston holding screw on the underside of the rod.
8. Click **Lower Syringe**.
9. After the syringe has been lowered, unscrew and remove the syringe from the valve.

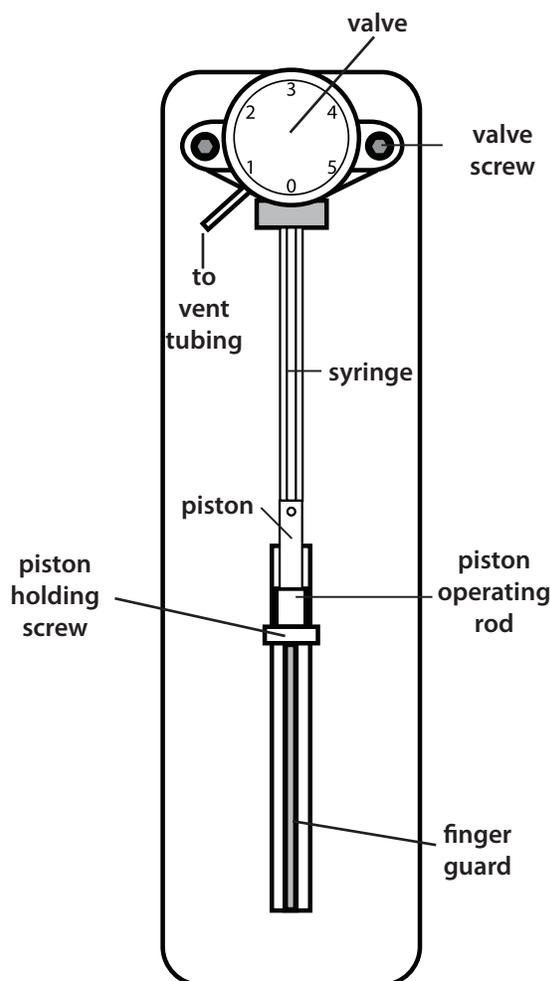


Figure 20: Syringe Installation

Clean the Syringe

After the syringe has been removed, it can be cleaned:

1. Place the syringe in a beaker containing methanol, and then aspirate and dispense several volumes of methanol through the syringe.
2. Place the syringe in a beaker containing distilled or deionized water, and then aspirate and dispense several volumes of water through the syringe.
3. Hold the syringe housing in one hand. Clean the syringe using a non-abrasive cloth dampened with alcohol. Remove the piston and clean the piston with a non-abrasive cloth dampened with alcohol.
4. Dry the syringe and piston using a clean, lint-free cloth.

Reinstall the Syringe

When the syringe is clean, reinstall it:

1. Fully tighten the syringe into the valve.
2. Click **Raise Syringe**.
3. Firmly tighten the piston holding screw to secure the syringe piston.

Valve

Clean the syringe pump valve with a nonabrasive cloth after any of the following situations have occurred:

- Corrosive or hazardous liquids have been pumped
- Possible back flow of liquids into the waste tubing
- Leakage
- System has not been used for a while

If the valve sticks, first try using the valve key (part number F123674) to turn the valve axle to possibly free the valve.

Remove the Valve

To clean the valve, first remove it from the syringe pump:

1. Disconnect all tubing from the valve.
2. Disconnect the syringe from the valve. Refer to the procedure for replacing the syringe on page 49.
3. Remove the two screws securing the valve to the syringe pump and then remove the valve.

Disassemble the Valve

1. Hold the valve body firmly in one hand. Using a 17 mm open-ended wrench, turn the valve axle guide counterclockwise and separate the two halves.
2. Pull the valve axle away from the valve main body.
3. Separate the ceramic stator from the ceramic rotor.

Note: Do not remove the ceramic stator from the valve main body.
4. Tap the valve axle guide against a solid level surface to remove the spring and PTFE end piece.

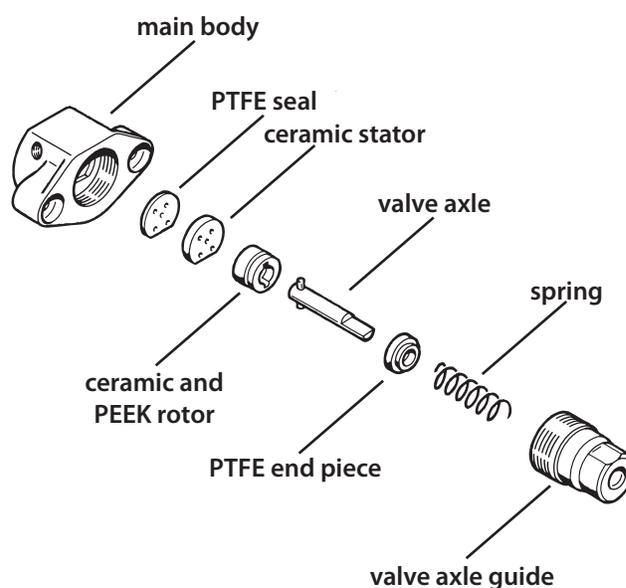


Figure 21: Valve (Disassembled)

Clean and Reinstall the Valve

1. Clean the disassembled parts of the valve using a non-abrasive cloth dampened with alcohol or by autoclaving.
2. Dry the components using a clean, lint-free cloth.
3. Reassemble the valve parts by reversing the above procedure.
4. Fully tighten the valve screws to secure the valve on the syringe pump.
5. Reinstall the syringe. Refer to the procedure for installing a new syringe on page 53.
6. Reconnect all tubing.

Part Replacement

Refer to the instructions in this section to replace the:

- **Syringe** on page 52
- **Valve** on page 53
- **Tubing** on page 52

Tubing

It is important to keep all tubing clean and free of crimps. Tubing that has become dirty, blocked, or crimped can result in poor accuracy and precision, or loss of air gap.

Replace both the transfer tubing and inlet tubing as needed. Refer to the **Parts and Accessories** appendix starting on page 61 for part numbers for replacement tubing. For tubing installation procedures, refer to **Plumbing Connections** starting on page 17.

Syringe

If necessary, refer to the diagram at right while performing the procedures below.

Remove the Syringe

The following procedures use TRILUTION® LH software.

1. Open a method with a VERITY 4060 Syringe Pump or VERITY 4260 Syringe Pump in the configuration.
2. Click **Run**.
3. Add a method to the sample list.
4. Select the Manual Control tab.
5. Select a method from the drop-down list of methods and then click **Go**. The instruments will initialize.
6. If using a VERITY® 4260 Dual Syringe Pump, select the syringe(s) that will be removed (Left Syringe and/or Right Syringe).
7. Disconnect the syringe piston from the piston operating rod by unscrewing the piston holding screw on the underside of the rod.
8. Click **Lower Syringe**.
9. After the syringe has been lowered, unscrew and remove the syringe from the valve.

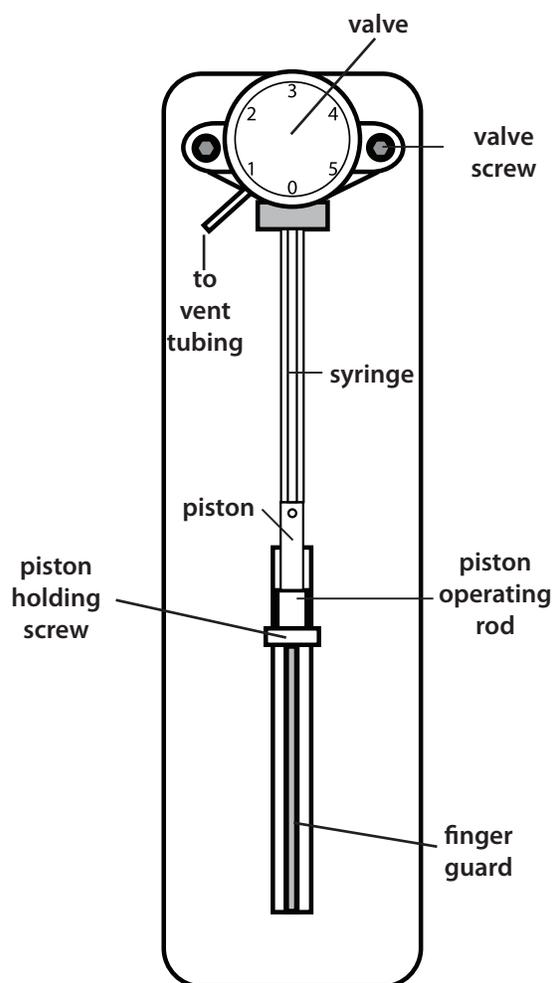


Figure 22: Syringe Installation

Install the New Syringe

1. Fully tighten the syringe into the valve.
2. Click **Raise Syringe**.
3. Firmly tighten the piston holding screw to secure the syringe piston.

Valve

To replace a valve on the syringe pump, follow the instructions below. If necessary, refer to the syringe pump diagram at right.

1. Disconnect the inlet, transfer, and vent tubing from the valve.
2. Disconnect the syringe from the syringe pump. Refer to the procedure for replacing the syringe on page 52.
3. Remove the two screws securing the valve to the syringe pump and then remove the valve.
4. Position the replacement valve and then fully tighten the valve screws to secure it.
5. Reinstall the syringe. Refer to the procedure for installing a new syringe on page 53.
6. Reconnect all tubing.

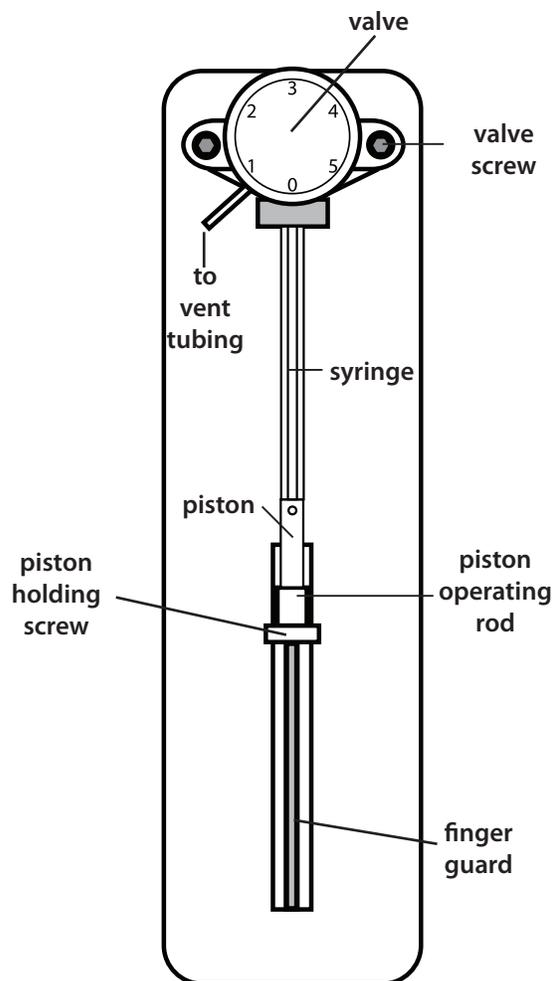


Figure 23: Syringe Installation

Troubleshooting

Chapter Five

This chapter provides information on the following topics:

- **Mechanical** on page 56
- **Electrical** on page 56
- **Hydraulic** on page 56
- **Error Codes and Messages** on page 58
- **Repair and Return Policies** on page 59

Mechanical

<i>Problem</i>	<i>Solution</i>
Valve does not fit on the syringe pump	Valve axle not properly aligned with the motor drive; ensure that the flat on the axle is aligned with the drive slot on the motor
Valve moves during aspirate or dispense	Valve mounting screw not tight; check that the mounting screws, syringe, and syringe mounting screw are all tight
Valve sticks	Clean the valve (refer to the instructions on page 51)

Electrical

<i>Problem</i>	<i>Solution</i>
Syringe pump does not respond	<p>Make sure power is turned on</p> <p>Check cabling connections</p> <p>Try a different AC outlet</p> <p>Restart the PC</p>

Hydraulic

<i>Problem</i>	<i>Solution</i>
Liquid leakage	Make certain all fittings are tight
Liquid leaving by the vent	<p>If tubing is kinked or blocked, replace tubing</p> <p>Ensure that the inlet tubing is clear</p> <p>Dirt may be trapped between the ceramic surfaces; clean the valve (instructions on page 51)</p>
Instrument will not draw in solvent	<p>Make certain all fittings are tight</p> <p>Check valve fitting threads on the syringe pump and replace if damaged</p>
No fluid being dispensed	<p>Make sure the syringe is tight within the valve fitting</p> <p>If tubing is kinked or blocked, replace defective tubing</p> <p>Replace the syringe pump valve if damaged</p>
Air gap breaks up	<p>When aspirating a liquid, if the air gap breaks up, check to see if the tubing is the correct size</p> <p>Reduce aspirate flow rate</p> <p>Increase volume of air gap</p> <p>Clean or replace any dirty tubing</p>

Problem	Solution
Syringe bubbles	<p>Make sure that all fittings are tight and air-free</p> <p>Make sure the syringe is tightened into the valve</p> <p>Clean the syringe if dirty (instructions on page 49)</p> <p>If any of the valve fittings are damaged, replace the valve</p>
Fluid leak	<p>Allow solvents to warm to room temperature before using</p> <p>Clean valve (instructions on page 51)</p> <p>Replace the syringe piston seal</p>
Incorrect aspirating and dispensing	<p>Check for leaks on all fittings</p> <p>Tighten or replace fittings on inlet and transfer tubing as needed</p> <p>Replace the valve if damaged</p> <p>Clean or replace transfer tubing</p>
Syringe stalls	<p>If the syringe stalls, inspect the tubing and valve for a blockage</p> <p>If the syringe stalls due to an accelerated aspirate or dispense rate, reduce the rate in the software</p>
Poor accuracy	<p>Clean or replace any dirty tubing</p> <p>Replace the syringe piston seal</p> <p>if the aspirate and dispense speeds are too fast, slow down the speeds to adapt to the tubing and probe type</p>

Error Codes and Messages

To obtain the error code and message, view the log file produced during the run in TRILUTION® LH:

To display the log file produced during a run:

1. Access the Run Results by clicking **Liquid Handling | Utilities | Run Results** or by clicking **Results** in the Application window.
2. Locate the run for which you want to view the log.
3. Click **View Log** or right-click on the Run and then select **View Log**. The file appears in a text editor box.

Refer to the table below for a list of the error codes and messages.

<i>Error</i>	<i>Error Text</i>	<i>Error</i>	<i>Error Text</i>
0	No Error	26	Invalid Valve Position
10	Unknown Command	27	Invalid Syringe Pump Volume
11	Invalid NV-RAM Address	28	Invalid Valve Move Speed
12	Emergency Stop Activated	29	Invalid Syringe Pump Flow Rate
13	Bad Parameter Entered	30	Invalid Valve Type
16	Character Limit	31	Invalid Syringe Size
18	Valve Park Location	32	Invalid Valve Selection
19	Syringe Pump Park Location	33	Invalid Syringe Pump Selection
20	Valve Unhomed	34	Missing Valve Encoder
21	Syringe Pump Unhomed	35	Missing Syringe Pump Encoder
22	Valve Moving	36	Set Pressure Offset to Zero
23	Syringe Pump Moving	37	Other Syringe Module in Error
24	Valve Stall	88	Error Unknown
25	Syringe Pump Stall		

Repair and Return Policies

Refer to the following information and then contact your local Gilson representative. Specific contact information can be found at www.gilson.com.

Before Calling Us

Your local Gilson representative will be able to serve you more efficiently if you have the following information:

- serial number and model number of the instruments involved
 - The serial number is located on the right side of the syringe pump.
- installation procedure you used
- list of concise symptoms
- list of operating procedures and conditions you were using when the problem arose
- list of all instruments in the configuration and the connections to those instruments
- list of other electrical connections in the room

Warranty Repair

Units covered under warranty will be repaired and returned to you at no charge. If you have any questions about applicability, contact your local Gilson representative.

Non-Warranty Repair

For out-of-warranty repairs, contact your local Gilson representative who will discuss service options with you and can assist in making arrangements to return the equipment, if necessary.

Return Procedure

Contact your local Gilson representative to obtain authorization before returning any Gilson equipment. To return a piece of equipment:

- Carefully pack the unit to prevent damage in transit. Check with your local Gilson representative regarding proper method of shipment. No responsibility is assumed by Gilson or your local Gilson representative for damage caused by improperly packaged instruments. Indicate the authorization on the carton and on the packing slip.
- Always insure for the replacement value of the unit.
- Include a description of symptoms, your name, address, phone number, and purchase order to cover repair costs, return and shipping charges, if your institution requires it.

Unit End-of-Life

When a unit reaches the end of its useful life, refer to www.gilson.com for directions and information on the end-of-life policy. This is in accordance with the European Union Directive 2002/96/EC on Waste Electrical and Electronic Equipment (WEEE).



Parts and Accessories

Appendix A

GX-Series ASPEC Systems

GX-241 ASPEC™ with VERITY® 4060 Single Syringe Pump

<i>Part Number</i>	<i>Description</i>
26150006	GX-241 ASPEC™ with VERITY® 4060 Single Syringe Pump

GX-271 ASPEC™ with VERITY® 4060 Single Syringe Pump

<i>Part Number</i>	<i>Description</i>
2614007	GX-271 ASPEC™ with VERITY® 4060 Single Syringe Pump

GX-271 ASPEC™ with VERITY® 4260 Dual Syringe Pump

<i>Part Number</i>	<i>Description</i>
2614008	GX-271 ASPEC™ with VERITY® 4260 Dual Syringe Pump

GX-274 ASPEC™ with Two VERITY® 4260 Dual Syringe Pumps

<i>Part Number</i>	<i>Description</i>
2614010	GX-274 ASPEC™ with two VERITY® 4260 Dual Syringe Pumps

Syringes

<i>Part Number</i>	<i>Description</i>
25025342	250 µL Syringe
25025347	500 µL Syringe
25025343	1 mL Syringe
25025344	5 mL Syringe
25025345	10 mL Syringe
25025346	25 mL Syringe

Piston Seals

<i>Part Number</i>	<i>Description</i>
F4015063	Piston Seal for 250 µL Syringe (pkg of 5)
F4015064	Piston Seal for 500 µL Syringe (pkg of 5)
F4015065	Piston Seal for 1 mL Syringe (pkg of 5)
F4015066	Piston Seal for 5 mL Syringe
F4015067	Piston Seal for 10 mL Syringe
F4015068	Piston Seal for 25 mL Syringe

Valve and Valve Parts

<i>Part Number</i>	<i>Description</i>
F4015040	Syringe Valve Assembly
F4015044	Rotor
F4015045	Seal
F4015046	Stator
F123674	Valve Key
49041019	Plug for Unused Ports
2502534811	Stainless Steel Thumbscrew for Syringe Piston

Tubing

Inlet

<i>Part Number</i>	<i>Description</i>
499484021	Inlet Tubing Assembly

Vent

<i>Part Number</i>	<i>Description</i>
F4420577	Vent Tubing

Plumbing Packages

<i>Part Number</i>	<i>Description</i>
2644700	Plumbing Package, GX-241/GX-271 ASPEC, 5 mL
2644701	Plumbing Package, GX-241/GX-271 ASPEC, 10 mL
2644702	Plumbing Package, GX-241/GX-271 ASPEC, 25 mL
2644703	GX-241/GX-271 ASPEC Air/Gas Plumbing Package
2644708	Dual Adapt Kit for Plumbing a GX-271 ASPEC™ with VERITY® 4260 Dual Syringe Pump
2644709	GX-271 ASPEC with Dual Syringe Pump Air/Gas Plumbing Package
2644704	Plumbing Package, GX-274 ASPEC, 5 mL
2644705	Plumbing Package, GX-274 ASPEC, 10 mL
2644706	Plumbing Package, GX-274 ASPEC, 25 mL
2644707	GX-274 ASPEC Air/Gas Plumbing Package

SPE Gas Pressure Regulator

<i>Part Number</i>	<i>Description</i>
25051376	SPE Gas Pressure Regulator Assembly 0–30 psi

Cables and Power Cords

<i>Part Number</i>	<i>Description</i>
7080316106	Power Cord, 220V
7080318107	Power Cord, 110V
32000012	USB Cable

Materials

Appendix B

Liquid Contact Materials

The information provided in the following table is accurate to the best of our knowledge and belief, but is intended for general information only.

<i>Material</i>	<i>Description</i>
Ekonol	<p>Ekonol has excellent solvent resistance with the exception of concentrated sulfuric acid and strong alkalis. The water absorption rate is low at 0.4% after 500 hours at 212° F. Ekonol Polyester is self-lubricating and provides excellent friction and wears properties.</p> <p>Ekonol Polyester is a very thermally stable polymer, making it easy to blend/fabricate with other high temperature materials. When combined with polytetrafluoroethylene (i.e., PTFE); it produces a composite material that has excellent temperature and wear resistance properties. The Ekonol Polyester/PTFE blend will not wear metal surfaces and resists self-wear better than any other PTFE composition. Applications for Ekonol Polyester/PTFE blends are varied and include packing sets, compressor ring sets, "O" ring seals, spring-loaded seals, lip seals, self-lubricating bearings and rotors or vanes of process pumps. Ekonol Polyester/PTFE works best under environmentally tough conditions where wear resistance, dimensional stability and corrosion resistance are critical.</p>
FEP	<p>Fluorinated ethylene propylene is another member of the fluorocarbon family with similar chemical properties. It is generally more rigid than PTFE, with somewhat increased tensile strength. It is typically more transparent than PTFE, slightly less porous, and less permeable to oxygen. FEP is not as subject to compressive creep at room temperature as PTFE, and because of its slightly higher coefficient of friction is easier to retain in a compression fitting.</p>

Liquid Contact Materials (Continued on page 66)

PEEK	<p>Considered relatively inert and biocompatible, polyetheretherketone tubing can withstand temperatures up to 100°C. Under the right circumstances, 0.005"–.020" ID tubing can be used up to 5000 psi for a limited time, and 0.030" to 3000 psi. Larger IDs are typically good to 500 psi. These limits will be substantially reduced at elevated temperatures and in contact with some solvents or acids.</p> <p>Its mechanical properties allow PEEK to be used instead of stainless in many situations and in some environments where stainless would be too reactive. However, PEEK can be somewhat absorptive of solvents and analytes, notably methylene chloride, DMSO, THF, and high concentrations of sulfuric and nitric acid. This tubing is highly prone to "kinking," or sealing off, if held in a sharp bend over time.</p>
PTFE	<p>Polytetrafluoroethylene is the generic name for the class of materials such as Teflon®. It offers superior chemical resistance but is limited in pressure and temperature capabilities. Because it's so easy to handle, it is often used in low pressure situations where stainless steel might cause adsorption. PTFE tubing is relatively porous, and compounds of low molecular weight can diffuse through the tubing wall.</p>

FEP, PEEK, and PTFE descriptions provided by Valco Instruments Company Inc (www.vici.com).

Ekonal description provided by Saint-Gobain Coating Solutions (www.coatingsolutions.saint-gobain.com)