

HOW DO I USE CONNECTED TOOLS AT THE BENCH?



1 	2 	3 	4 	5 
CHECK the activity and calibration status of your device in PipetteScope®	REQUEST a pipette calibration/service if needed	OPEN the PipettePilot® app	LOAD your protocol from SciNote or your tablet	EDIT and save changes to your protocol
6 	7 	8 	9 	10 
CONNECT your PIPETMAN® M Connected and environment sensor to the app	PLACE your plate support and microplate on the screen	FOLLOW the light on the app that shows where, when and how to pipette	REVIEW your automatically generated report and add comments	SAVE the report locally or send it to your ELN